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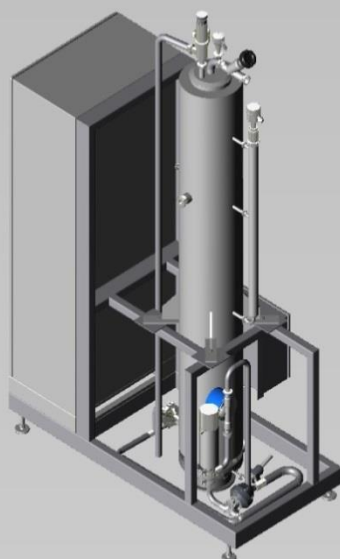
Pure Steam Generation

PureSteam

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The Principle

Pure steam is used for sterilization of tanks, pipe-works and filling machines, but also for direct injection into a product. For pure steam generation the feed water enters a tubular heat exchanger with an expansion chamber above. In the tubular heat exchanger the pre-treated water is heated and steam is generated. The tubes of the heat exchanger are connected with the expansion chamber that serves as a separating column. For production of pyrogen-free steam the chamber contains a demistor. After entering the expansion chamber, the steam moves further upwards in the column. By gravity and due to the specific column design, unevaporated water droplets fall back into the lower section of the chamber, thus separating from the steam. Since the droplets may contain contaminants like endotoxins, it is vital that they are not carried forward with the steam. The conductivity of the pure steam condensate can be monitored. Optionally, the unit is expandable to a distillation system for combined generation of pure steam and pure water.

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The Centec Group

Centec offer fully automated, skid-mounted process units and high precision measurement technology from a single source. Our systems and sensors are engineered to perfectly meet the most demanding requirements of the brewery, beverage, food and pharmaceutical industries. With a team of experienced engineers, we aim to create quantifiable added value for our customers. We are your partner from planning and design through to commissioning of your plant.

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